SmartScope iGO Quick Start Guide

iTunes Install:

- 1. Visit the iTunes Store and search for the APP "Wi-Viewer"
- 2. Install "Wi-Viewer"

Android Install:

- 1. Visit the Google Play Market and search for the APP "Mi-Cam"
- 2. Install "Mi-Cam"
- A. Open the battery door on the SmartScope iGO and make note of the SSID.
- B. Turn the SmartScope iGO on and leave on for about 30 seconds.
- C. Turn the tablet on and go to SETTINGS and choose WiFi.
- D. Choose the network SSID from the inside battery cover.
- E. Password is "12345678"
- F. After connected, go to the HOME screen and open the application.

***** iOS devices ONLY: Touch the ON button on the bottom left of the screen to get the live image.



Snapshot Button to take a picture.

Begin with focus dial turned all the way to the right. Slowly turn to the left until the object comes into focus.

If touching the object with the lens tip, continue to turn the focus dial to the left and the image will come into focus a second time for even greater magnification.

ON / OFF Switch

Light Intensity Dial



*****Technical support, please call 281-312-1297, 1-866-722-2051 or email support@smartschoolsystems.com

